

Priyanka Sharma

M.Sc Final
(Chemistry)


Organic Chemistry

Index

Chemistry

Sr. No.	Name of the Experiment	Page No.	Date of Experiment	Date of Submission	Remarks
1.	To prepare benzilic acid from benzoin				<u>Suven</u>
2.	To prepare benzanilide from benzophenone by the backmann-rearrangement.				
3.	To prepare 2-phenyl indole from phenyl hyrazine by Fisher-indole synthesis				
4.	To separate & identify the given mix. containing one liquid & two solid comp. & pre. their suitable derivatives.				
5.	To separate & identify given mix. containing one liquid & two solids comp. & prepare their suitable derivatives.				<u>Suven</u>
6.	To separate & identify the given mixture containing one liquid & two solid comp. prepare their suitable derivatives				<u>Suven</u>

Chemical Rxn -

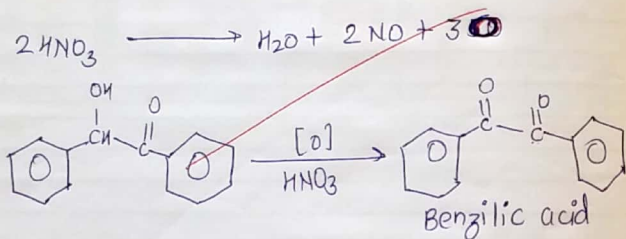

Date _____

Expt. No. _____

Page No. _____

Experiment = 1

Object:

To prepare benzilic acid from benzoin

Step I → Preparation of benzil from benzoin

Chemical Requirement -

- Benzoin → 8.5 gm
- Glacial acetic acid → 40ml
- Conc. HNO_3 → 20ml
- Ice


Procedure -

- In a round bottom flask place 8.5 gm of the benzoin, 40ml of glacial acetic acid & 20ml conc. HNO_3 .
- Heat the mixture on a steam bath for two hours. Cool the flask in an ice bath add 150ml of water, mix thoroughly & allow the yellow ppt of benzil to settle.
- Filter the product with solution, wash thoroughly with water & dry the product by pressing the crystal with a clean glass stopper.
- Although again the benzil obtained is sufficiently pure for conversion to derivatives or to benzilic acid, it may be purified by crystallisation from methanol.

A.P.P.P.

Teacher's Signature _____

Chemical Rxn

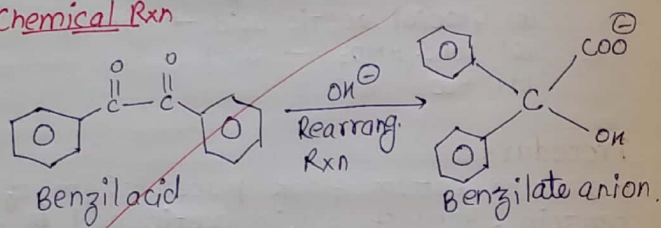

Expt. No. _____

Date _____

Page No. _____

Result -

Appearance \rightarrow Yellow solid Crystal
Yield \rightarrow 8 gm
M.P \rightarrow 95°C.

Step - II \rightarrow Preparation of benzilic acid from benzil.

Chemical Requirement -

- (i) Benzil from I-step
- (ii) Potassium hydroxide - 5ml
- (iii) Ethyl alcohol - 15 ml
- (iv) Conc. HCl
- (v) gce.

Procedure -

- (i) Dissolve 5 gm of Potassium hydroxide pellets in 10 ml of water in a 100 ml flask.
- (ii) Add 15 ml of ethyl alcohol & mix well the resulting solution.
- (iii) Add 5 gm of pure benzil to the resulting solution.
- (iv) Fit a reflux water condenser to the flask & boil the contents on a water bath for 10-15 minutes.
- (v) Transfer the contents of the flask to a small beaker or a porcelain dish, cover with a watch glass all it to stand undisturbed for several hours preferably overnight until potassium benzilate is completely crystallised.

ASAP

Teacher's Signature _____

Filtered the product & wash it with ice cold ethanol.

Dissolve the potassium benzilate in about 50ml of water & then dissolve or add concⁿ solⁿ of HCl drop by drop till the precipitation of the free acid is complete.

Filter the product wash with water press dry & recrystallise from boiling water or benzene.

Result:


Appearance — Colourless shining crystal

Yield — 4.5 gm.

M.P — 150°C.

Silviana
28/4/18

Chemical Rxn -

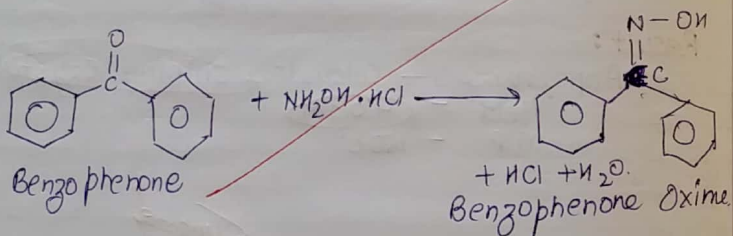

Expt. No. _____

Date _____

Page No. _____

Experiment - 2

Object:

To prepare benzanilide from benzophenone by the ~~backmann-rearrangement~~.

~~Step-I: Preparation of benzophenone oxime from benzophenone~~

Chemical Requirement -

- (i) Benzophenone - 5 gm
- (ii) Hydroxylamine hydrochloride - 3 gm
- (iii) Rectified spirit - 10 ml
- (iv) Sodium hydroxide (Pelets) - 5.6 gm
- (v) Conc. HCl - 15 ml


Procedure -

- (i) Place 5 gm of benzophenone, 3 gm of hydroxylamine hydrochloride, 10 ml of rectified spirit & 2 ml of water in a flask.
- (ii) Now add 5.6 gm of NaOH portion wise & with shaking & cool the flask. In a case than Rxn became to vigorous.
- (iii) Boil the solⁿ under reflux condenser for 5 min.
- (iv) Cool the solⁿ pour the contents to a solⁿ of 15 ml of conc. HCl in 100 ml of water in a 500 ml beaker.

APAR

Teacher's Signature _____

Chemical Rxn. :-

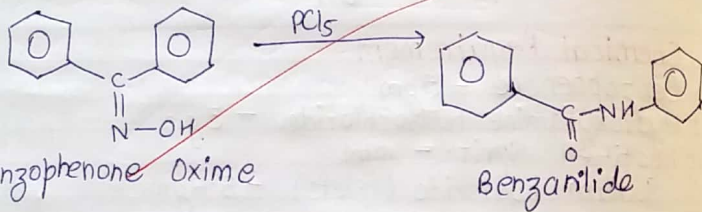

Result-

Appearance \rightarrow Crystalline Solid
Yield - 5 gm
M.P - 142°C

Step-II: Preparation of benzalidine from Benzophenone oxime by backmann's Rearrangement.

Chemical properties-

- (i) Benzophenone Oxime - 4 gm
- (ii) Potassium penta chloride - 6 gm.
- (iii) Anhydrous Ether - 40 ml


Procedure:

- (i) Dissolve 4 gm of benzophenone oxime in 40 ml of anhydrous ether in a conical flask.
- (ii) Add 6 gm of powdered PCl_5 & 6 ml of pure thionyl chloride to the Rxn mixture.
- (iii) Shake well the contents & distil off the solvent on water bath.
- (iv) Cool the residue, add 50 ml water & boil the resulting solⁿ for several min.

Result:

Appearance = White Crystalline Solid
Yield = 3 gm
M.P = 169°C

Chemical Rxn -

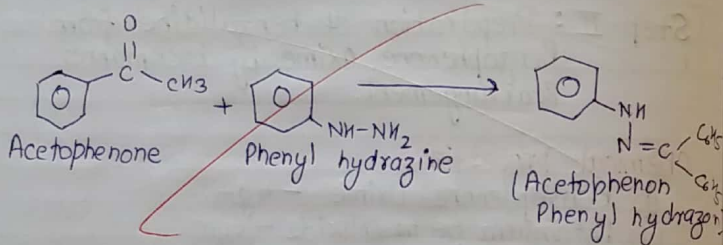

Date _____

Expt. No. _____

Page No. _____

Experiment - 3

Object-

To prepare 2-Phenylindole from phenyl
hydrazine by Fisher-Indole synthesis

Step-I - Preparation of Phenylhydrazin
derivative of acetophenone.

Chemical required-

- (i) Acetophenone - 10ml
- (ii) Phenyl hydrazin - 12ml
- (iii) Ethanal - 30ml
- (iv) Acetic acid - 2-3ml

Procedure-

- (i) In a round bottom flask. Place 10ml of acetophenone, 12ml Phenyl hydrazin, 30 ml ethanal containing 2-3 drops of glac. CH_3COOH .
- (ii) The solⁿ was warmed at 100°C for 15 min
- (iii) After the poured the mix. into a beaker containing ice cool water.
- (iv) The phenyl hydrazine derivative was filtered & washed with dil HCl.

Result: Brown colour crystal form acetophenone
phenyl hydrazone
Yield = 7.3 gm M.P = 106°C

Teacher's Signature _____

Step - II : To prepare 2-Phenyl indole from acetophenone Phenyl hydrazine.

Chemical requirement -


- (i) Acetophenone phenyl hydrazine - 3gm
- (ii) Polyphosphoric acid - 40ml
- (iii) Water - 100ml.

Procedure -

Acetophenone phenyl hydrazone formed in step - I was heated with Polyphosphoric acid in boiling oil with 15min in round bottom flask.

Result :

Crystal of 2-Phenyl indole
Yield = 4.89 gm
M.P = 187°C

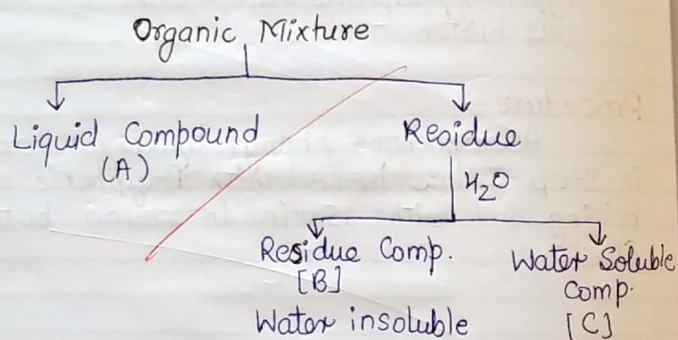

Expt. No. _____

Date _____

Page No. _____

Experiment - 4 (1)

Object:

To separate & identify the given mix. containing one liquid & two solid compound & prepare their suitable derivative.

Separation:

The given solid-liquid mixture are separated by distillation at 80°C liquid compound. A solid mix. of concⁿ B & C was separated by H_2O .

[A] Identification of Compound "A":

Primary test:

- (i) Physical State - liquid
- (ii) Colour - Colourless
- (iii) Odour - Specific odor
- (iv) Solubility - Insoluble in water


Observation Table:

S.No.	Experiment	Observation	Inference
①	Ignition test: Compound take in spectula	Burn without smoky flem	Aliphatic
②	Litmus test: Litmus paper dip in compound sol ⁿ	No effect on litmus paper	Compound is neutral.

APR 2019

Teacher's Signature _____

Derivative of acetone.

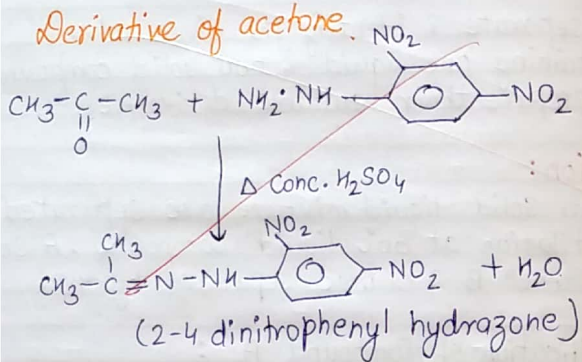

③ Element test:

(a) N-Test: L.S + few drops (1-2) of (NaOH Solⁿ + freshly prepared FeSO₄ Solⁿ + Δ + cool + add conc. H₂SO₄) Green/blue ppt is not obtained. N-absent

(b) S-Test:

L.S + dil CH₃COOH + Few drops of (CH₃COOH)₂ Pd. Black ppt. is not obtained. S-absent

(c) X-test:

L.S + dil HNO₃ + AgNO₃ Solⁿ White & Yellow ppt is formed. X-halogen is present

④ Functional gp. test

Compⁿ + 2ml alcohol + (1-2) Pellets KOH + Δ + cool + dil HNO₃ + AgNO₃ Solⁿ White & Yellow ppt. is formed. X-halogen is present

5 Determination of boiling point

B.P of Compound 61°C Chloroform may be

6 Compound test:

(i) Comp. + Crystal of resorcinol + aq. NaOH + Δ Red colour is obtained. Chloroform Confirms

JKK

Teacher's Signature _____

B7 Identification of Solid Comp. (B) :


Primary test:

- (i) Physical State - Solid Crystalline
- (ii) Colour - Yellow.
- (iii) Odour - Specific Colour
- (iv) Solubility - Insoluble in water

Observation Table -

S.No	Experiment	Observation	Inference
1	Ignition test: Comp ⁿ take on spectula	Compound burns with smoky flame	Comp ⁿ is aromatic
2	Litmus test: Comp ⁿ Sol ⁿ + Litmus paper	No effect on the litmus paper	Comp. is neutral.
3	Element test:		
(a)	N-test : L.S + NaOH. Sol ⁿ + FeSO ₄ Sol ⁿ + Δ + Cool + Conc. H ₂ SO ₄	Green & blue ppt. is not obtained	N is absent.
(b)	S-test : L.S + dil CH ₃ COOH + (CH ₃ COO) ₂ Pd Sol ⁿ	Black ppt. is not obtained	"S" is absent
(c)	X-test : L.S + conc. HNO ₃ + AgNO ₃ Sol ⁿ	White ppt is not obtained.	X-absent.

Derivative :->

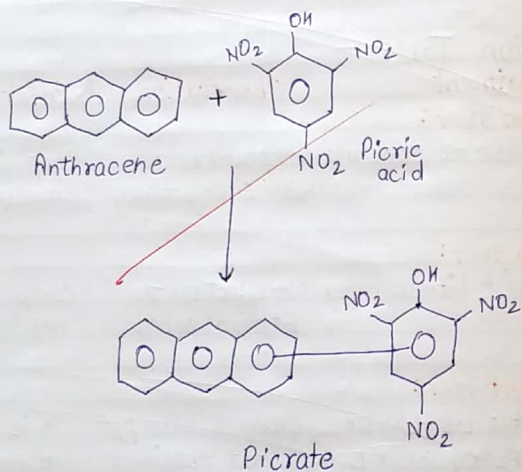

Expt. No. _____

Date _____

Page No. _____

4. Functional gp. test:

Substrate + benzen solⁿ + dil solⁿ of iodine + shake

Solⁿ remains violet

Aromatic hydrocarbon confirms

5. Melting pt.

M.P of the compound

216°C

Anthracene may be

6. Compound test:

Compⁿ + Substrate solⁿ of picric acid in the benzene + Δ

Deep red colour of solid is obtain

Anthracene is confirmed.

7. Derivative -

Compⁿ + ethanol + saturated solⁿ of Picric acid + Δ + cool + Filter

Red colour solid is obtained

Picrate

8. M.P of derivative

M.P of Picrate

138°C

Picrate confirmed

(c) Identification of Solid Compound 'c' :-

Primary test:

- (i) Physical state - Crystalline Solid
- (ii) Colour - Brown
- (iii) Odour - Carbolic Odour
- (iv) Solubility - Poorly soluble in water

ASAP

Teacher's Signature _____

(4) Functional gp. test-

(i) Compⁿ + FeCl₃ Solⁿ Violet colour obtained Phenol may be, +nt.

(ii) Libermann-nitrous test:


(a) Comp. + Na Nitrate + conc. H₂SO₄ Green colour Confirms.

(b) Dil it with H₂O Red colour obtain

(c) Then add it NaOH. It convert into blue colour Phenol Confirms.

(5) Determination of m.p.:

M.P of compound 110°C

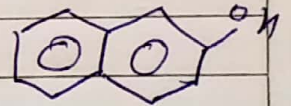

β -Naphthol may be.

(6) Compound test:


(i) Comp. + NaOH solⁿ + CCl₄ + Δ Blue colour is obtained β -naphthanol confirms.

Derivative -

Comp. + benzene + Picric acid + Shake + crystal is brown warm benzene.

Yellow crystal obtained

Picrate

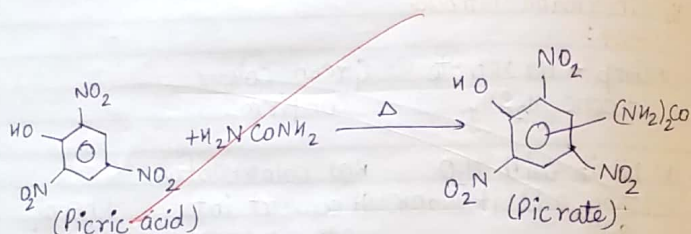

M.P of Picrate 156°C Picrate Confirmed

Result:

In the given organic mixture the following comp. are present.

Compound - A:


Comp. (A) is chloroform ($CHCl_3$)
 B.P of Comp - 61°C
 Derivative - 2,4 dinitro derivative

Compound - B:

Comp. (B) is anthracene (c1ccc2ccccc2c1)
 M.P of Comp - 216°C
 Derivative - Picrate
 M.P of derivative - 138°C

Compound - C:

Comp. "C" is β -naphthol (Oc1ccc2ccccc2c1)
 M.P of Comp. - 110°C
 Derivative - Picrate
 M.P of derivative - 156°C

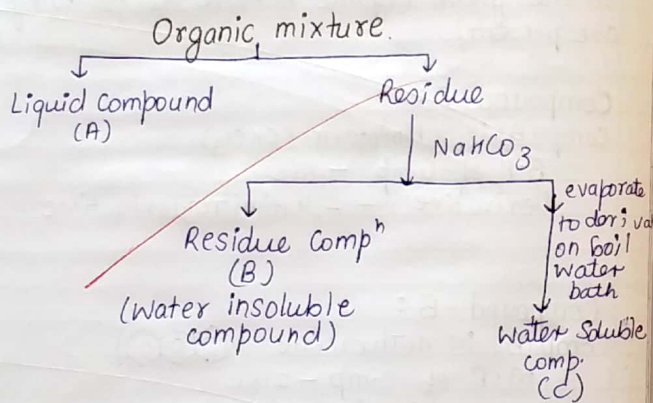

Expt. No. _____

Date _____

Page No. _____

Experiment = 5

Object:

To separate & identify the given mix. containing one liquid & two solid compⁿ & prepare their suitable derivatives.

Separation:

The given solid-liquid mix. are separated by distillation of 80°C liq. compound A & solid mix. of compⁿ B & C was separated by H₂O.

Identification of Compound (A):

Primary test:

- ① Physical state - liquid
- ② Colour - Colourless
- ③ Odour - Specific odour
- ④ Solubility - Soluble in water

Observation Table:

S.No	Experiment	Observation	Inference
	Comp. take in the Spetulla	Comp. ⁿ burn with smoky flame	Comp. ⁿ is aliphatic

Teacher's Signature _____

Teacher's Signature _____

Litmus Test:Litmus paper dip
in compⁿ solⁿNo effect on the
litmus paperComp. is
Neutral**Element test:**(a) **N-test:** L.S + NaOH +
FeSO₄ + Δ + cool + conc.
H₂SO₄Green & blue ppt/
Solⁿ not obtained

N-absent

(b) **S-test:**L.S + dil CH₃COOH +
(CH₃COO)₂ Pd solⁿBlack ppt is not
obtained

S-absent


(c) **X-test:**L.S + dil HNO₃ + AgNO₃ solⁿWhite ppt. is not
obtained

X-absent

Functional gp. test:-**Ceric ammonium mixture
test:**Comp. + Few drops ceric
ammonium nitrate +
ShakeRed colour
obtainedAlcohol gp
may be**Determination of B.P.:**B.P of compⁿ

78°C

Ethyl alcohol
may be

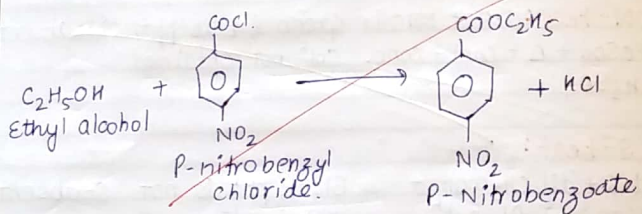

Expt. No. _____

Page No. _____

Compound test :

Comp. + I_2 Solⁿ + NaOH + Δ . Yellow colour obtained Ethyl alcohol confirm.

Derivatives -

Alcohol + p-nitro benzoyl Chloride + Δ + NaHCO₃ Solⁿ + Cool. Solid p-nitro benzoate is obtained P-nitrobenzoate.

M.P.

57°C

Identification of Solid Comp(B) -

Primary Test :

- ① Physical state - Crystalline solid
- ② Colour - White
- ③ Odour - specific odour
- ④ Solubility - Soluble in hot water

Observation Table -

S.No.	Experiment	Observation	Inference
①	Ignition test: Comp. take in the Spectulla	Comp. burns with smoky flame	Comp. is aromatic

Teacher's Signature _____

Teacher's Signature _____

Litmus test:

Comp. Sol ⁿ + litmus paper	No effect	Comp. is neutral
---------------------------------------	-----------	------------------

Element test:**(a) N-test:**

L.S + NaOH + FeSO ₄ Sol ⁿ + Δ + Cool + Conc. H ₂ SO ₄	Green/blue ppt. / Sol ⁿ are obtain	N - present
---	---	-------------

(b) S-test:

L.S + dil CH ₃ COOH + (CH ₃ COO) ₂ Pd Sol ⁿ	Black ppt is not obtained	S-absent
---	---------------------------	----------

(c) X-test:

L.S + conc. HNO ₃ + AgNO ₃ Sol ⁿ	White ppt. is not obtained	X-absent
---	----------------------------	----------

Functional gp. Test:

Comp. + H ₂ O + Shake + H ₂ O ₂ Sol ⁿ + Δ + Cool + FeCl ₃ Sol ⁿ	Red/blue colour obtain	Aromatic amide may be
---	------------------------	-----------------------

M.P of Compound
M.P of Compound

129°C

Benzamide may be

Compound Test:

Comp. + dil NaOH + Δ + cool + dil HCl	White ppt. is obtain	Benzamide is confirm
---------------------------------------	----------------------	----------------------

Derivations :

Comp. + aniline White ppt. obtain
+ Δ

M.P of derivative 162°C Benzaldehyde confirms,

Result:

In the organic mix. following comp. are +nt

Compound A :

Comp. (A) is ethyl alcohol

B.P of Comp - 78°C

Derivatives - p-nitrobenzoate

M.P of Derivatives - 57°C


Compound B :

Comp. (B) is Pthalic acid


M.P of the comp. - 195°C

Derivative - Amide

M.P of derivative - 220°C

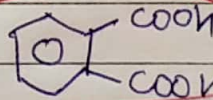**Compound C :**

Comp. (C) is benzamide

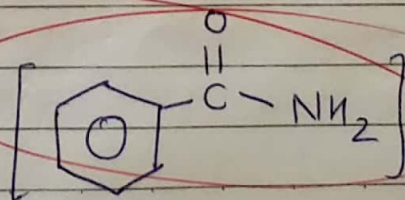


Date _____

Expt. No. _____

Page No. _____

M.P of Comp. -129°C
Derivative - Benzaldehyde
M.P of derivative -162°C .

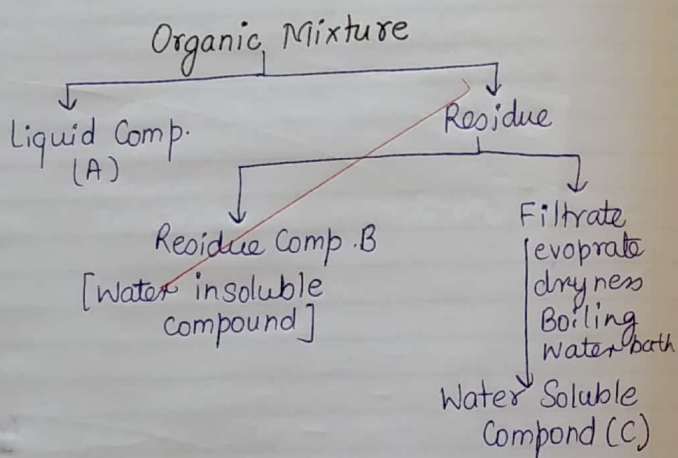

Expt. No. _____

Date _____

Page No. _____

Experiment = 6

Object:

To separate & identify the given mixture containing one liquid & two solid compⁿ prepare their suitable derivatives.

Separation:

The given solid-liquid mixture are separated by distillation at 80°C liquid comp [A] & solid mix. of comp. (B) & (C) was separated by H₂O.

Identification of Compound (A):

Primary test:

- ① Physical state - liquid
- ② Colour - Colourless
- ③ Odour - Specific odour
- ④ Solubility - Soluble in water

Observation table:

S.No.	Experiment	Observation	Inference
1)	Ignition test: Comp. take in the spectulla	Comp. burns with non-smoky flame	Comp. is aliphatic

Teacher's Signature _____

Teacher's Signature _____

Litmus test:Litmus paper dip
in comp. solⁿ

No effect

Comp. is
neutral**Element test:****(a) N-test:**L.S + few drops of
NaOH + freshly prep.
FeSO₄ + Δ + Cool +
Conc. H₂SO₄~~Green/blue ppt.~~
is not obtained

N-absent


(b) S-test:dil CH₃COOH + (CH₃COO)₂Pd
SolⁿBlack ppt.
not obtained.

S-absent

(c) X-test:L.S + dil HNO₃ + AgNO₃
SolⁿWhite ppt. not
obtainedX-absent
~~Alcohol may~~
be.**Functional gp. test:****Ceric ammonium
Nitrate**Comp. + Ceric ammo-
-nium nitrate + ShakeRed colour
obtainedalcohol
may be**Melting pt. of Compound:**M.P of compⁿ

65°C

methyl
alcohol
may be

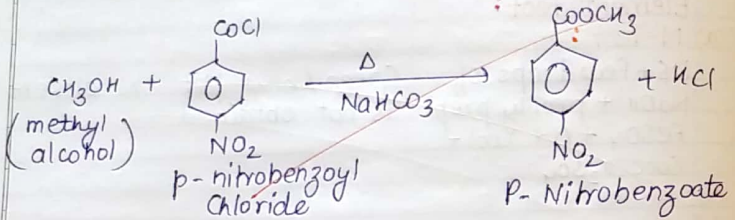

Expt. No. _____

 Date _____
 Page No. _____

Compound test:

Comp. + Salicylic acid + conc. $\text{H}_2\text{SO}_4 + \Delta + \text{H}_2\text{O}$	Smell of winter green oil	CH_3OH confirms
--	---------------------------	---------------------------------

Derivatives:

Alcohol + p-nitrobenzoyl Chloride + Δ + NaHCO_3 + Cool	Crystal of p-nitrobenzoate obtain	p-nitrobenzoate confirmed
--	-----------------------------------	---------------------------

Determination of M.P.:

M.P. of Comp.	96°C	p-nitrobenzoate confirmed
---------------	------	---------------------------

Identification of Solid Compound 'B': -
Primary test:

- ① Physical State - Crystalline solid
- ② Colour - Specific odour
- ③ Odour - Colourless
- ④ Solubility - Soluble in water

Observation table:

S.No.	Experiment	Observation	Inference
	Ignition test: Comp. taken in the spectulla	Burns with non-smoky flames	Comp. is aliphatic

Teacher's Signature _____

P.P.P.

Litmus test:Litmus paper dip
in comp. SolⁿBlue change in
redacidic in
nature**Element test:****(a) N-test:**L.S + few drops of
+ NaOH Solⁿ + FeSO₄
Solⁿ + Δ + Cool +
conc. H₂SO₄~~Blue / Green ppt.~~
is not obtain

N-absent

(b) S-test:L.S + CH₃COOH +
(CH₃COO)₂Pd SolⁿBlack ppt. is
not obtain

S-absent

(c) X-test:L.S + dil. HNO₃ +
AgNO₃ SolⁿWhite ppt. is
not obtain

X-absent

(4) Functional Gp. test:**= Sodium bicarbonate****test:**Comp. + NaHCO₃
+ Δ CO₂ gas evolved-COOH gp.
may be.**(5) M.P of Compound:**

M.P of Comp.

185°C

Succinic
acid may
be

Compound test:

Comp. + Resorcinol + conc. H_2SO_4 + Δ ,	Red-brown colour obtain.	
--	-----------------------------	--

To mix + NaOH,	Orange-green colour obtain	Succinic acid confirmed
----------------	-------------------------------	----------------------------

Derivatives :

Comp. + $SOCl_2$ + Δ + Cool + conc. NH_3 Sol ⁿ + cool it once cube	White crystal obtain	Amide confirm
--	-------------------------	------------------

Determination of M.P.:

M.P of derivative comp.	260°C	Amide confirms
----------------------------	-------	-------------------

Identification of Solid Compound (C) :**Primary test:**

- ① Physical State - Solid
- ② Colour - Colourless
- ③ Odour - Specific odour
- ④ Solubility - Insoluble in water

Observation Table :

S.No.	Experiment	Observation	Inference
	Ignition test: Comp. take in spectula	Comp. burns with smoky flames	Comp. is aromatic

Litmus Test:Litmus paper dip
in comp Solⁿ

No effect

Comp. is
neutral in nature**Element test:****(a) N-test:**L's + few drops of
NaOH Solⁿ + freshly
prepared FeSO₄ Solⁿ
+ Δ + cool +
Conc. H₂SO₄Green/blue
ppt. Solⁿ is
obtained

N-present

(b) S-test:L's + CH₃COOH +
(CH₃COO)₂ Pd SolⁿBlack ppt is
not obtain

S-absent

(c) X-test:L's + dil. HNO₃ +
AgNO₃ SolⁿWhite ppt. is
not obtain

X-absent

Functional gp. test:**Mulliken & Baker
test:**Alcohol + CaCl₂ + Zinc
dust powder + Δ +
cool + tollen reagentBrown/black
colour ppt.
ObtainNitro gp.
may be**M.P of compound.**

M.P

90°C

m-dinitroben-
zene may
be

Compound test:

(i) Comp. + NaOH + Δ	Voilet colour	m-dinitro-
(ii) Comp. + dil NaOH + Δ glucose	obtain	-benzene confirms

Determination of M.P:

M.P of derivative comp.	90°C	m-dinitro-benzene confirms
-------------------------	------	----------------------------

Result:

In the following compound (organic mix.) present -

Compound (A) -

Compound (A) is methanol (CH_3OH)

M.P - 65°C

Derivative - p-nitrobenzoate

M.P of derivative - 96°C

Compound (B) -

Compound (B) is succinic acid


M.P - 185°C

Derivative - Amide

M.P of derivative - 260°C

Compound - "C":

Comp. "C" is dinitrobenzene

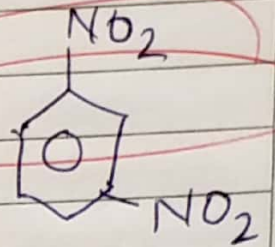

M.P of Comp. -90°C .

~~Shreya~~
28/4/18